



SPME Applications

Amines, Aromatics, Explosives

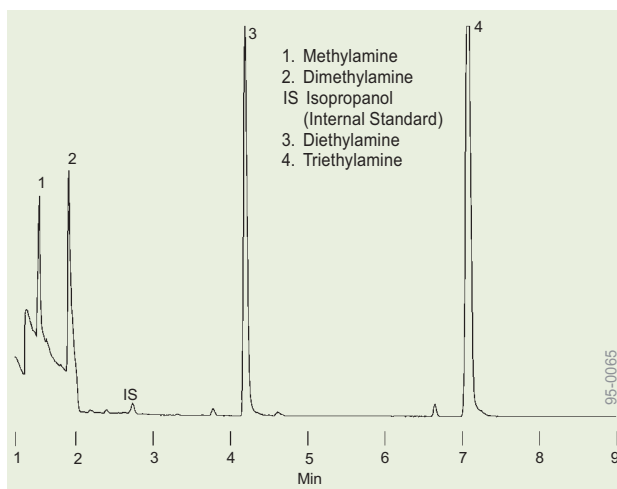

Amines (SPME/GC)


Sample: amines at 5ppm in water, 27% NaCl, pH 9.5
 SPME Fiber: polydimethylsiloxane/divinylbenzene, 65µm film
 Cat. No.: 57310
 Extraction: 20 min, fiber immersed in water, rapid stirring
 Desorption: 5 min, 270°C
 Column: SPB-1 SULFUR, 30m x 0.32mm ID, 4.0µm film
 Cat. No. 24158
 Oven: 50°C (2min) to 150°C at 10°C/min
 Carrier: helium, 25cm/sec (set at 50°C)
 Det.: GC/MS, selected ions at 0.6 sec/scan
 Inj.: splitless/split, closed initial 5 min, 270°C, 0.75mm ID splitless liner

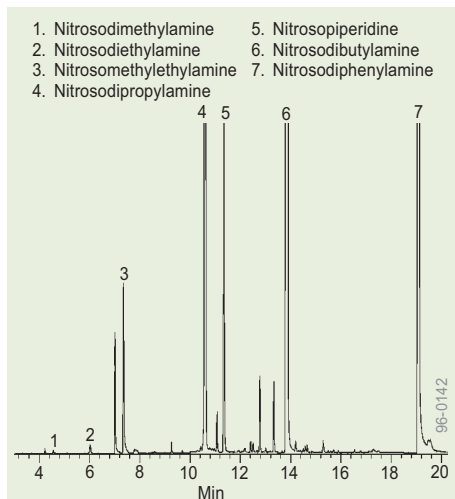

Chromatogram courtesy of J. Clark, Liggett Group, Inc.

Nitrosamines (SPME/GC)

Sample: analytes in water + 25% KCl, pH 10.0
 SPME Fiber: polydimethylsiloxane/divinylbenzene, 65µm film
 Cat. No.: 57310-U
 Extraction: immersion, 15min (rapid stirring)
 Desorption: 270°C, 1min
 Column: PTA-5, 30m x 0.32mm ID, 0.5µm film
 Cat. No.: 24331
 Oven: 50°C (1min) to 250°C at 10°C/min, hold 2 min
 Carrier: helium, 30cm/sec
 Det.: GC/MS (quadrupole, selected ion mode)
 Inj.: splitless, 250°C (0.75mm ID liner)

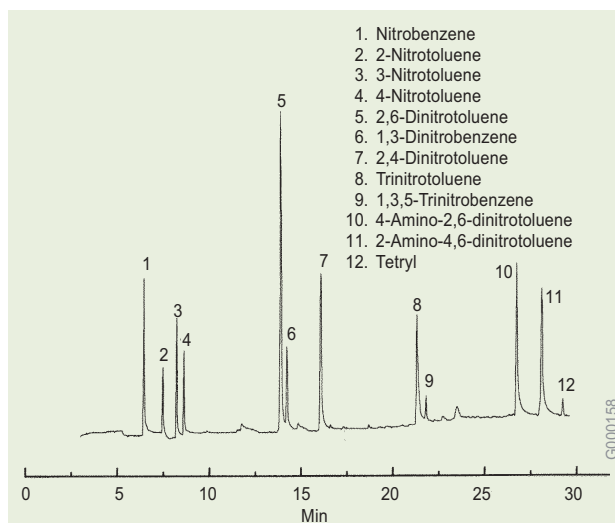

Reproduced from The Analysis of High Explosives Residues by Solid-Phase Microextraction Followed by HPLC, GC/ECD and GC/MS by R. Almirall, Metro-Dade Police Crime Laboratory, Miami, FL, US. Presented at Academy of Forensic Science 2/13/98.

Explosives in Water (SPME/GC)

Sample: 50ng/mL each explosive in water
 SPME Fiber: polydimethylsiloxane/divinylbenzene, 65µm film
 Cat. No.: 57310-U
 Sampling: immersion, 30 min
 Desorption: 5 min, 250°C
 GC Column: cyanopropyl silicone, 30m x 0.25mm ID, 0.25µm film
 Supelco
 Equivalent: SPB-1701 (Cat. No. 24113)
 Oven: 95°C (3 min) to 182°C at 8°C/min (4 min) to 250°C at 8°C/min (6 min)
 Carrier: nitrogen, 60mL/min
 Det.: ECD, 250°C
 Inj.: split/splitless, 180°C


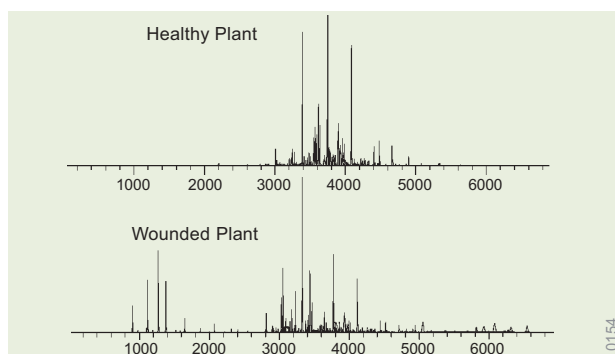


Figure provided by A.-K. Borg-Karlson, The Royal Institute of Technology, Department of Chemistry, Stockholm, Sweden.

Mono- and Sesquiterpenes in Spruce Seedlings (SPME/GC)

Sample: headspace surrounding *Picea abies* seedlings
 SPME Fiber: polydimethylsiloxane, 100µm film
 Cat. No.: 57300-U
 Extraction: headspace, 60 min
 Desorption: 30 sec, 225°C
 Column: poly(5% diphenyl/95% dimethylsiloxane), 30m x 0.25mm ID, 0.25µm film
 Supelco
 Equivalent: Equity-5 (Cat. No. 28089-U)
 Oven: 40°C (4 min) to 200°C at 4°C/min, hold 16 min
 Carrier: helium, 10psi
 Det.: EID, 70ev
 Inj.: splitless (splitter opened after 30 sec), 225°C

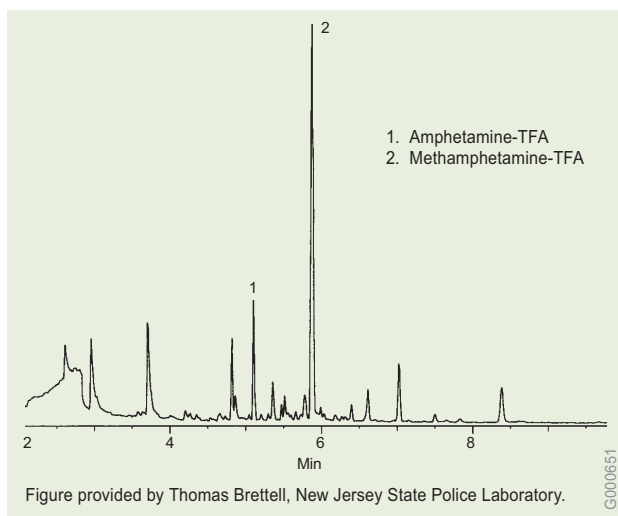

Amphetamines (SPME/GC)

Sample: 1mL urine + 0.7g K₂CO₃ in 20mL headspace vial, equilibrated at 80°C, 30 min

SPME Fiber: polydimethylsiloxane, 100µm film
Cat. No.: 57300-U

Extraction: headspace, 3-5 min, 80°C

Derivatization: methyl bis-trifluoroacetamide (headspace, 0.5 min, ambient)

Desorption: 1 min, 270°C


Column: methylsiloxane, 12.5m x 0.2mm ID, .033µm film

Supelco
Equivalent: Equity-1 (available on request)

Oven: 60°C (1 min) to 140°C (4 min) at 30°C/min, then to 276°C at 20°C/min, 4 min

Det.: MS, full scan

Inj.: splitless (closed 1 min), 270°C

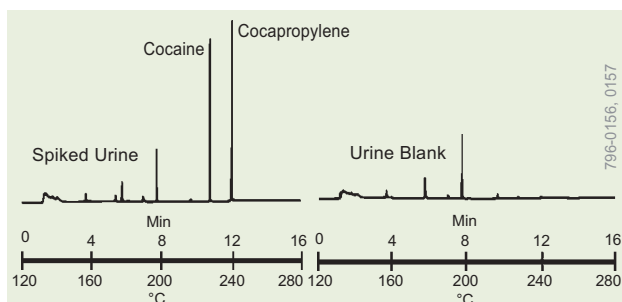

Cocaine in Urine (SPME/GC)

Sample: 0.5mL urine (250ng each analyte, 20µL 2.5% NaF) in 1mL vial

SPME Fiber: polydimethylsiloxane, 100µm film
Cat. No.: 57300-U

Extraction: immersion, 30 min

Desorption: 3 min, 240°C

Column: polydimethylsiloxane, 30m x 0.32mm ID, 0.25µm film


Supelco
Equivalent: Equity-1 (Cat. No. 28055-U)

Oven: 120°C to 280°C at 10°C/min

Carrier: helium, 3mL/min

Det.: NPD, 280°C

Inj.: splitless (splitter opened after 1 min), 240°C

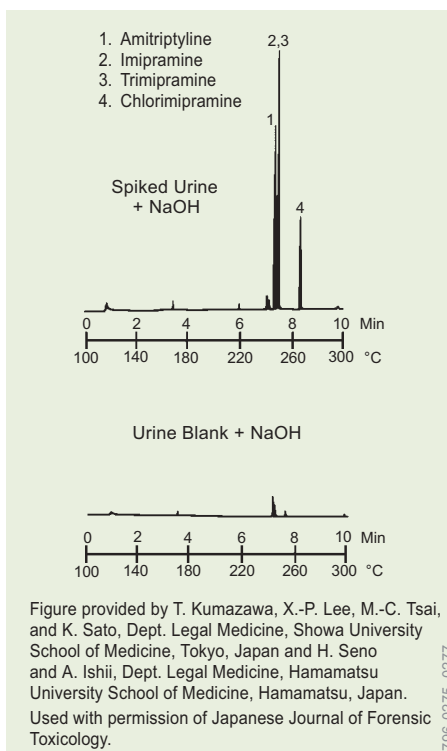

Tricyclic Antidepressants in Urine (SPME/GC)

Sample: 1mL urine (1µg each analyte + 50µL 5M NaOH) in 7.5mL vial

SPME Fiber: polydimethylsiloxane, 100µm film
Cat. No.: 57300-U

Extraction: headspace, 15 min, 100°C (sample incubated 30 min)

Desorption: 3 min, 280°C

Column: polydimethylsiloxane, 30m x 0.32mm ID, 0.25µm film

Supelco
Equivalent: Equity-1 (Cat. No. 28055-U)

Oven: 100°C to 300°C at 20°C/min


Carrier: helium, 3cm/sec

Det.: FID, 280°C

Inj.: splitless (splitter opened after 1 min), 280°C


SPME Applications

Foods, Flavors and Fragrances

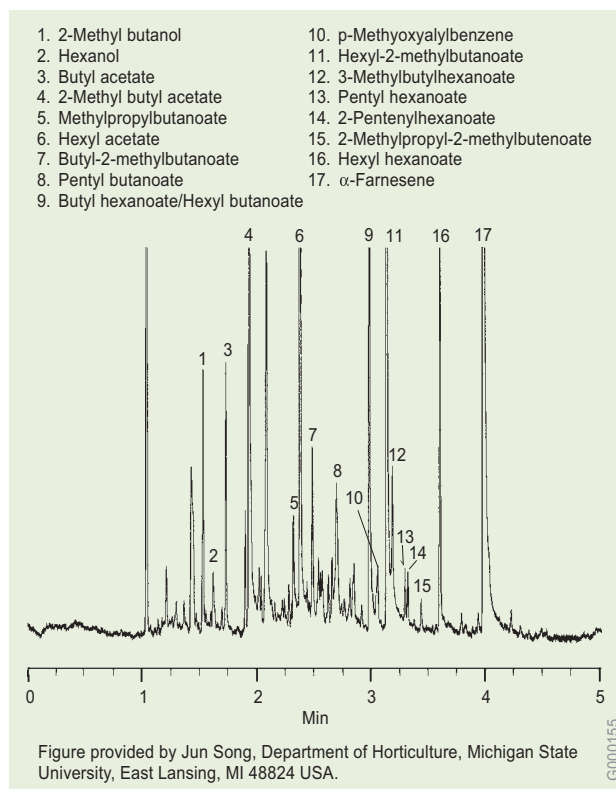

Aroma Volatiles from Stored Apples (SPME/GC)

- Sample: 300-450g intact fruit in a 3 liter flask at ambient temp., exposed to an air flow of 25-30mL/min.
- SPME Fiber: polydimethylsiloxane/divinylbenzene, 65 μ m film headspace extraction (as described for Sample), 4 min; analytes desorbed (250°C) to a cool (<100°C) precolumn, then volatilized by heating to 150°C in 15 sec.
- Cat. No.: 57310-U
- GC Column: poly(5% diphenyl/95% dimethylsiloxane) type, 25m x 0.25mm ID, 0.34 μ m film
- Supelco Equivalent: Equity-5 (available on request)
- Oven: 40°C (1.5 min) to 250°C at 50°C/min, hold 1 min
- Carrier: helium, 1.5mL/min (44cm/min)
- Det.: mass spectrometer, m/z = 40-300 (40 spectra/sec)

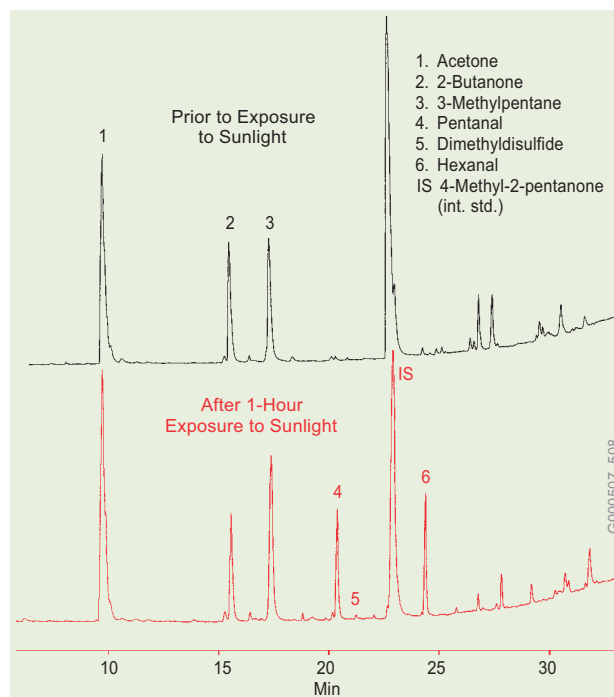

Milk Sample Off-Flavors (SPME/GC)

- Sample: 3g of 2% milk + 10 μ L internal standard solution (20 μ g/mL 4-methyl-2-pentanone) (9mL GC vial)
- SPME Fiber: PDMS/Carboxen, 75 μ m film
- Cat. No.: 57318
- Extraction: headspace, 15 min with constant stirring at 45°C
- Desorption: 5 min, 250°C
- Column: Supel-Q PLOT, 30m x 0.32mm ID
- Cat. No.: 24242
- Oven: 70°C (2 min) to 140°C at 6°C/min (2 min hold) then to 220°C at 6°C/min (5 min hold)
- Carrier: helium, 35cm/sec
- Det.: GC/MS ion trap, m/z = 33-300

Acid

- | | |
|----------------|--------------------|
| 1. Acetic | 7. Hexanoic |
| 2. Propionic | IS 2-Ethylhexanoic |
| 3. Isobutyric | 8. Heptanoic |
| 4. Butyric | 9. Octanoic |
| 5. Isovaleric | 10. Nonanoic |
| 6. Valeric | 11. Decanoic |
| 12. Undecanoic | 13. Dodecanoic |

Free Fatty Acids in Parmesan Cheese (SPME/GC)


- Sample: 100mg cheese in 40mL vial
- SPME Fiber: 65 μ m Carbowax/divinylbenzene StableFlex
- Cat. No.: 57336-U
- Sampling: 65°C 15 min (headspace)
- Desorption: 250°C, 1 min
- GC Column: Nukol, 15m x 0.25mm, 0.25 μ m film
- Cat. No.: 24106-U
- Oven: 50°C (2 min) to 220°C at 10°C/min
- Carrier: helium, 30cm/sec
- Det.: FID, 260°C
- Inj.: splitless/split (closed 1 min), 250°C

SPME Applications Foods, Flavors and Fragrances

Coffee Components*

- | | | |
|----------------------------|--------------------------------|------------------------------------|
| 1. 2-Methyl furan | 17. Methyl pyrazine | 31. 2-Ethylpyrazine |
| 2. 2-Butanon | 18. 4-Methyl thiazole | 32. 2-Ethyl-3,5-dimethylpyrazine |
| 3. 2-Pentanone | 19. 3-Hydroxy butanon | 33. Glycerol |
| 4. 3-Methyl butanal | 20. Dimethyl phenol (isomer) | 34. 2,3-Diethylpyrazine |
| 5. 2,5-Dimethylfuran | 21. 1,2-Ethanediol monoacetate | 35. 2-Ethyl-3,6-dimethylpyrazine |
| 6. 2-Acetyloxy-2-propanone | 22. 2,5-Dimethylpyrazine | 36. 2-Furancarboxaldehyde |
| 7. 2-Ethyl hexanol | 23. 2,3-Dimethylpyrazine | 37. 2-Isopropenylpyrazine |
| 8. Dimethylsulfoxide | 24. 2-Ethylpyrazine | 38. 3,5-Diethyl-2-methylpyrazine |
| 9. Phenol | 25. 2,6-Dimethylpyrazine | 39. Furfural formate |
| 10. Hexanal | 26. 2-Ethyl-6-methylpyrazine | 40. 2-Furonyl ethanone |
| 11. 2-Methyl thiophene | 27. 2-Ethyl-5-methylpyrazine | 41. Methyl benzoylformate |
| 12. n-Methyl pyrrole | 28. Trimethylpyrazine | 42. Furanmethanol acetate |
| 13. 4-Methylphenol | 29. 2-Ethyl-3-methylpyrazine | 43. 5-Methyl-2-furancarboxaldehyde |
| 14. 2-Ethyl pyrrole | 30. 2,6-Diethylpyrazine | 44. Furanmethanol propionate |
| 15. Pyridine | | |
| 16. Pyrazine | | |

- | |
|--|
| 45. Furfanyl furan |
| 46. Pyridine methanol |
| 47. 2-Methyl-5-propenylpyrazine |
| 48. Furanmethanol |
| 49. 3-Ethyl-4-methyl-2,5-furandione |
| 50. Pyrazine carboxamide |
| 51. 2-Ethyl-3-hydroxy-4H-pyran-4-one |
| 52. 1-(2-Furanylmethyl)pyrrole |
| 53. 2-Methoxyphenol |
| 54. 1-(1H-Pyrrole-2-yl)ethanone |
| 55. 4-Ethyl-2-methoxy phenol |
| 56. 3-Phenylpropenal or 2-Methylbenzofuran |
| 57. 3,5-Dimethylbenzoic acid |

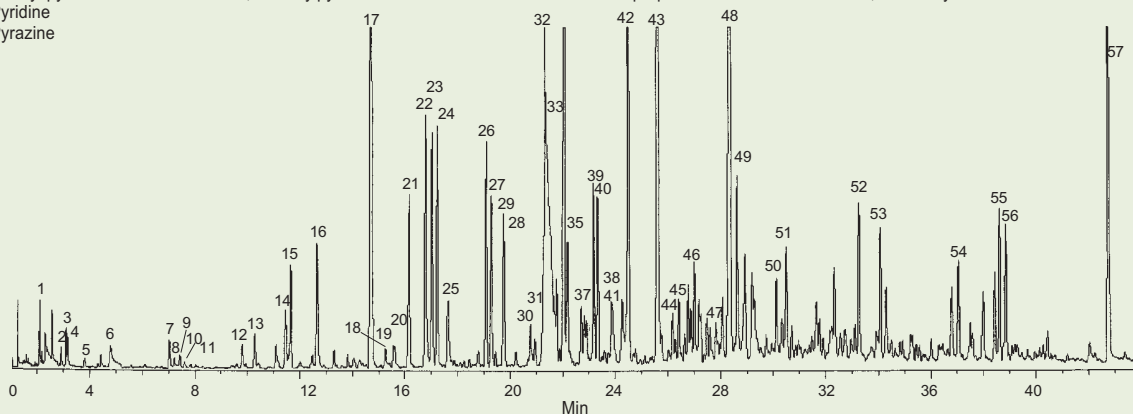

* Identification of some of the pyrazines is not absolute.

Ground Coffee (SPME/GC)

Sample: 5g ground coffee in 40mL vial
 SPME Fiber: 50/30µm DVB/Carboxen/PDMS (StableFlex fiber)
 Cat. No.: 57348-U
 Extraction: 30 min, 65°C (headspace)
 Desorption: 270°C, 5 min

GC Column: SUPELCOWAX 10, 30m x 0.25mm ID, 0.25µm film
 Cat. No.: 24079
 Oven: 40°C (5 min) to 230°C at 4°C/min
 Det.: ion trap MS, m/z = 30-350, 0.6 sec/scan
 Inj.: splitless/split (closed 0.5 min), 270°C (0.75mm ID inlet liner)


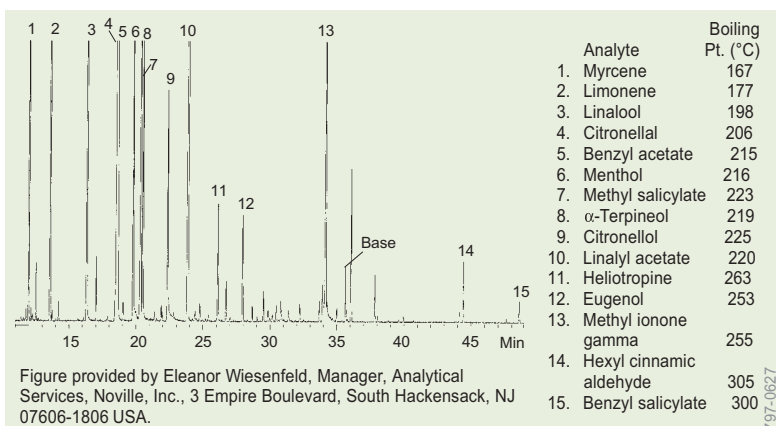


Figure provided by Eleanor Wiesenfeld, Manager, Analytical Services, Noville, Inc., 3 Empire Boulevard, South Hackensack, NJ 07606-1806 USA.

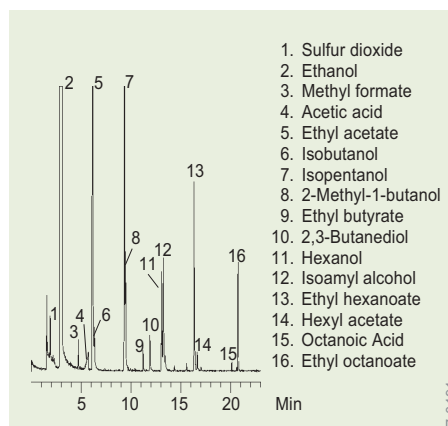

Volatiles in White Wine (SPME/GC)

Sample: White wine + 25% NaCl
 SPME Fiber: Carboxen/PDMS
 Cat. No.: 57318
 Extraction: 10 min, 40°C (headspace)
 Desorption: 3 min at 290°C
 Column: VOCOL, 30m x 0.25mm ID, 1.5µm film
 Cat. No.: 24205-U
 Det.: GC/MS, quadrupole, m/z = 31-240


Shampoo Base with Fragrance (SPME/GC)

Sample: 0.5g shampoo base with 1% fragrance test mix, diluted 1:1 with NaCl/water in a 10mL vial
 SPME Fiber: polydimethylsiloxane, 100µm film
 Cat. No.: 57300-U
 Extraction: headspace, 5 min, 80°C (with sample stirring)
 Desorption: 5 min, 250°C
 Column: methyl silicone, 50m x 0.25mm ID, 0.25µm film
 Supelco
 Equivalent: Equity-1 (available on request)
 Oven: 35°C to 240°C at 3°C/min
 Cryo: -50°C w/liquid CO₂
 Flow: 0.5mL/min
 Det.: MSD, scan m/z = 29-350 AMU
 Inj.: split, 250°C

SPME Applications

Foods, Flavors and Fragrances

- | | |
|--------------------|----------------------|
| 1. Acetic acid | 12. Heptanoic acid |
| 2. Pentanal | 13. Octanal |
| 3. Butanoic acid | 14. Octanoic acid |
| 4. Propyl acetate | 15. Nonanone |
| 5. Methyl butyrate | 16. Nonanal |
| 6. Hexanal | 17. Butyl hexanoate |
| 7. Octane | 18. Decanal |
| 8. Methyl hexanal | 19. Undecanone |
| 9. Hexanoic acid | 20. Pentyl hexanoate |
| 10. Heptanone | 21. Dodecanone |
| 11. Heptanal | 22. Methyl heptanol |
| | 23. Dodecanal |

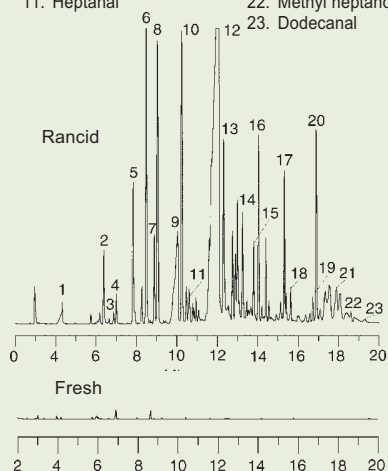

G000075,76

Potato Chips (SPME/GC)


Sample: 3g crushed potato chips in 15mL vial
 SPME Fiber: DVB/Carboxen/PDMS, 50/30µm film, StableFlex
 Cat. No.: 57328-U
 Extraction: headspace, 65°C, 20 min, with stirring
 Desorption.: 3 min, 250°C
 GC Column: SPB-1 SULFUR30m x 0.32mm ID, 4.0µm film
 Cat. No.: 24158
 Oven: 45°C (1.5 min) to 250°C at 12°C/min, hold 10 min
 Carrier: helium, 40cm/sec
 Det.: quadrupole mass spectrometer, m/z = 35-290 at 0.6 sec/scan

Some Volatile Components in Peanut Butter

- Carbon disulfide
- 3-Methylbutanal
- Pentanal
- Dimethyl disulfide
- Hexanal
- 1-Methyl pyrrole
- 4-Methyl-pentene-2-one
- Heptanal

Pyrazines in Peanut Butter

- 2-Methyl pyrazine
- 2,5-Dimethyl pyrazine
- 2,3-Dimethyl pyrazine
- 2-Ethyl pyrazine
- 2,6-Dimethyl pyrazine
- 2-Ethyl-6-methyl pyrazine
- 2-Ethyl-5-methyl pyrazine
- Trimethyl pyrazine
- 2-Ethyl-3-methyl pyrazine
- 2,6-Diethyl pyrazine
- 2-Ethyl-3,5-dimethyl pyrazine
- 2,3-Diethyl pyrazine
- 2-Methyl-5-isopropyl pyrazine
- 3-Ethyl-2,5-dimethyl pyrazine
- 5-Methyl-2-propyl pyrazine
- 2-Methyl-5-propyl pyrazine
- 2-Ethenyl-6-methyl pyrazine
- 3,5-Diethyl-2-methyl pyrazine
- 2-Ethenyl-5-methyl pyrazine
- 2-Methyl-6-cis propenyl pyrazine
- 2-Allyl-5-methyl pyrazine

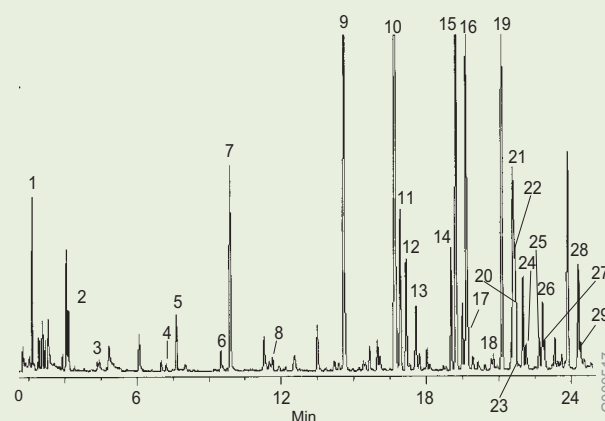

G000517

Pyrazines in Peanut Butter (SPME/GC)

Sample: 5g peanut butter in 40mL vial
 SPME Fiber: divinylbenzene-Carboxen-polydimethylbenzene StableFlex
 Cat. No.: 57328-U
 Extraction: headspace, 30 min at 65°C in heating block
 Desorption: 270°C for 5 min
 Column: SUPELLOWAX 10, 30m x 0.25mm, 0.25µm film
 Cat. No.: 24079
 Oven: 40°C (5 min) to 230°C at 4°C/min
 Inj.: splitless/split, closed 0.5 min, 270°C, with 0.75mm liner
 Det.: ion trap mass spectrometer, m/z = 30-350 at 0.6 sec/scan
 Selected ions used for quantitation.

SPME Applications Hydrocarbons and Pesticides


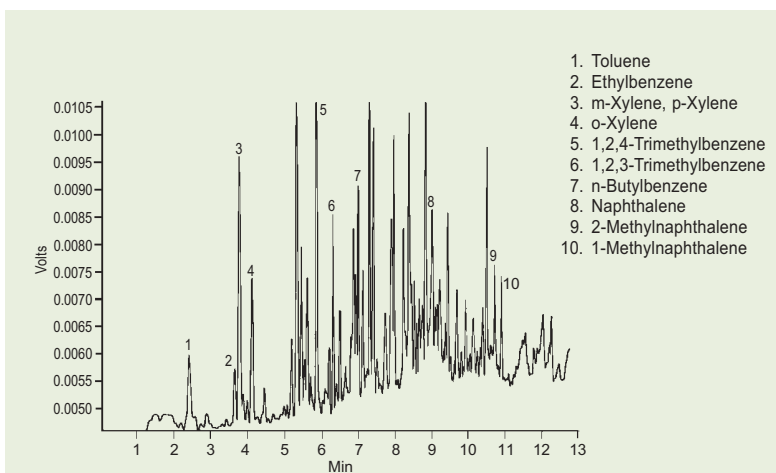


Figure courtesy of José Almirall, Crime Laboratory Bureau, Metro-Dade Police Department, Miami, FL, USA, and Kenneth Furton and Juan Bruna, Department of Chemistry, Florida International University, Miami.

Reproduced from the Journal of Forensic Sciences. Copyright American Society for Testing and Materials. Reprinted with permission.

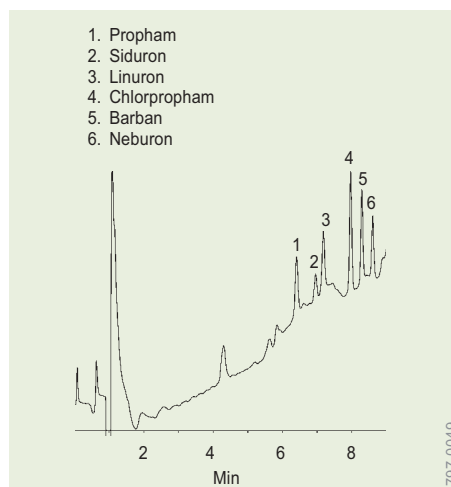

Carbamate and Urea Pesticides (SPME/HPLC)

Sample: 3mL water containing 8ng/mL of each analyte in 10% NaCl

SPME Fiber: PDMS/DVB, 60µm film

Cat. No.: 57317

Extraction: immersion, 40 min, rapid stirring

Desorption: static, 5 min in acetonitrile:water (65:35);
dynamic, valve open during run

Column: SUPELCOSIL LC-8, 15cm x 4.6mm ID,
5µm particles

Cat. No.: 58220-U

Mobile Phase: acetonitrile:water
(18:82 to 65:35 in 9 min, hold 3 min)

Flow Rate: 2.0mL/min

Temp.: 35°C

Det.: UV, 240nm

Gasoline in Arson Samples (SPME/GC)

SPME: polydimethylsiloxane phase fiber, 100µm film
headspace sampling (20 min) 10 sec desorption (splitless mode)

Cat. No.: 57300-U

Column: polydimethylsiloxane phase, 30m x 0.25mm ID, 0.25µm film

Supelco


Equivalent: Equity-1 (Cat. No. 28046)

Oven: 35°C (2 min) to 220°C at 10°C/min, hold 2 min,
to 300°C at 30°C/min, hold 5 min

Carrier: helium, 1mL/min (split 50:1)

Det.: FID, 300°C

Inj.: splitless (closed 3 min), 220°C (2mm ID injector liner)

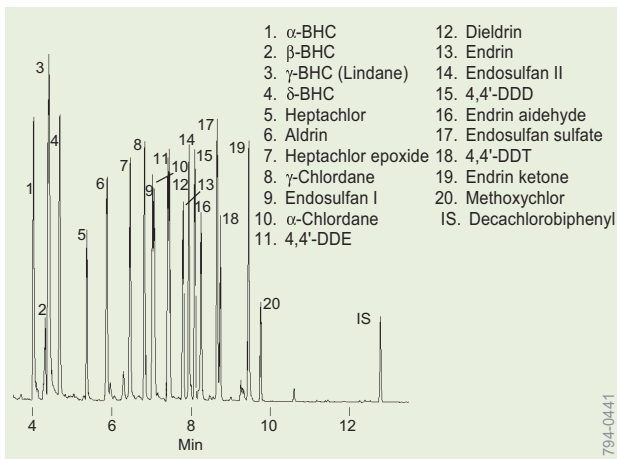

Chlorinated Pesticides (SPME/GC)

Sample: 200ppt each in 2mL water

Extraction: 15 min immersed in water with rapid stirring

SPME Fiber: polydimethylsiloxane, 100µm film

Cat. No.: 57300-U

Column: SPB-5, 15m x 0.20mm ID, 0.20µm film


Cat. No.: 24165-U

Oven: 120°C (1 min) to 180°C at 30°C/min, then to 290°C at 10°C/min

Carrier: helium, 37cm/sec (set at 120°C)

Det.: ECD, 300°C

Inj.: 260°C (splitless - closed 3min)

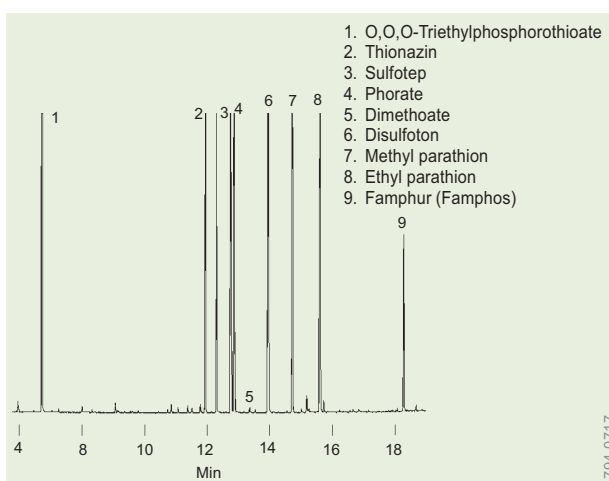

Organophosphorous Pesticides (SPME/GC)

Sample: 50ppb each in 1.5mL saturated salt water, pH 7.2

Extraction: 20 min immersion with rapid stirring

SPME Fiber: polydimethylsiloxane, 100µm film

Cat. No.: 57300-U

Column: PTE-5, 30m x 0.25mm ID, 0.25µm film

Cat. No.: 24135-U

Oven: 60°C (1 min) to 300°C at 12°C/min, hold 5 min


Carrier: helium, 30cm/sec

Det.: MS (m/z = 45-400, 0.6sec/scan)

Inj.: splitless (0.75mm ID liner), 270°C


SPME Applications

Herbicides, PCBs, and Phenols

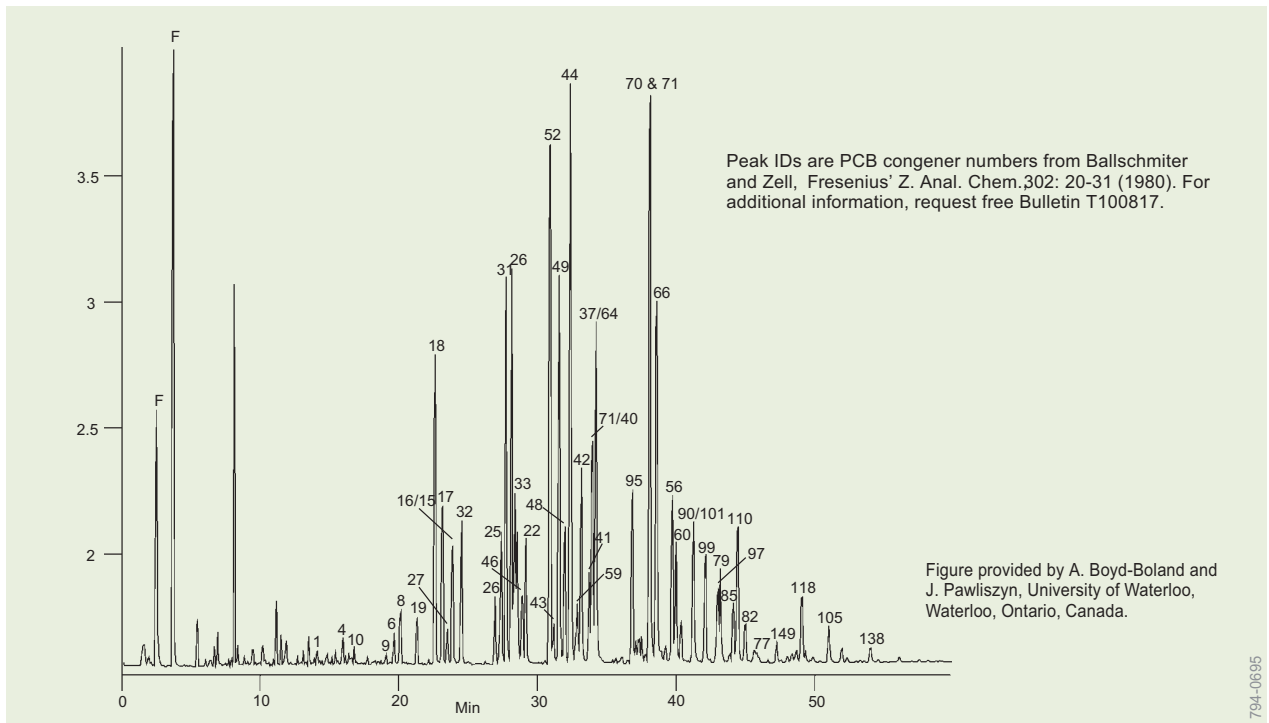

PCB Congeners in Stream Sediment (SPME/GC)


SPME Fiber: PDMS, 100 μ m film
 Cat. No.: 57300-U
 Extraction: headspace, 90°C/30 min
 Desorption: 280°C/2 min, splitless
 Column: SPB-Octyl, 30m x 0.25mm ID, 0.25 μ m film
 Cat. No.: 24218-U
 Oven: 50°C/1 min at 10°C/min to 150°C/0 min
 Sample: 5g stream sediment, downstream from industrial sites

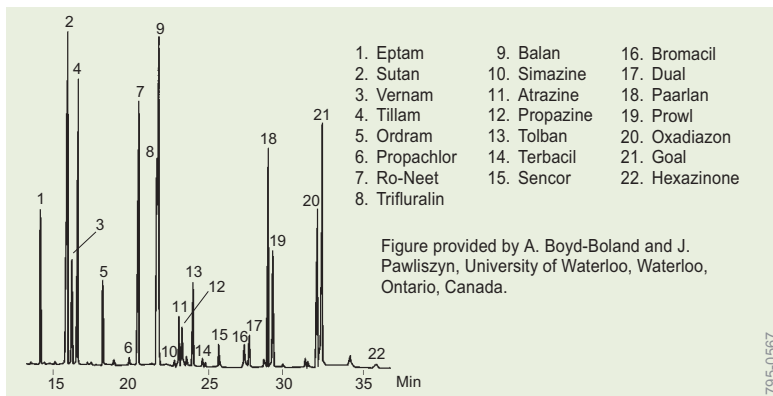

Nitrogen-Containing Herbicides Extracted from Water (SPME/GC)

Sample: 4mL (water + 100ng/mL each analyte 1g/mL NaCl), pH 2, in 4.6mL vial
 SPME Fiber: polyacrylate, 85 μ m film
 Cat. No.: 57304
 Extraction: immersion, ambient temp., 50 min (constant stirring)
 Desorption: 5 min, 230°C
 Column: PTE-5, 30m x 0.25mm ID, 0.25 μ m film
 Cat. No.: 24135-U
 Oven: 40°C (5 min) to 100°C at 30°C/min, to 275°C at 5°C/min
 Carrier: helium, 40cm/sec, set at 40°C
 Det.: MS (Ion Trap Manifold: 250°C;
 Mass Scan Range: m/z = 45-400 at 0.6 sec/scan)
 Inj.: splitless, 230°C

IS 2-Fluorophenol (int. std.)	10. 4-Chloro-3-methylphenol (int. std.)
1. Phenol	11. 2,4,5-Trichlorophenol
2. 2-Chlorophenol	12. 2,4,6-Trichlorophenol
3. 2-Methylphenol	13. 2,4-Dinitrophenol
4. 3-Methylphenol	14. 4-Nitrophenol
5. 4-Methylphenol	15. 2,3,4,6-Tetrachlorophenol
6. 2-Nitrophenol	16. 2-Methyl-4,6-dinitrophenol
7. 2,4-Dimethylphenol	IS 2,4,6-Tribromophenol (int. std.)
8. 2,4-Dichlorophenol	17. Pentachlorophenol
9. 2,6-Dichlorophenol	18. Dinoseb

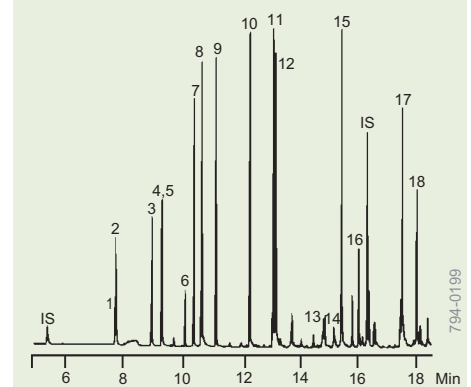


Phenols in Water (SPME/GC)

Sample: 50ppb phenols in 1.8mL saturated salt water, pH 2
 Extraction: 20 min immersed in water with rapid stirring
 SPME Fiber: polyacrylate, 85 μ m film
 Cat. No.: 57304
 Desorption: splitless, 280°C (closed 3 min)
 Column: PTE-5, 30m x 0.25mm ID, 0.25 μ m film
 Cat. No.: 24135-U
 Oven: 40°C (4 min) to 260°C at 12°C/min
 Carrier: helium, 40cm/sec at 40°C
 Det.: MS, Scan Range m/z = 45-465 at 0.6 sec/scan


SPME Applications

Polymers and Oligomers, Polynuclear Aromatic Hydrocarbons

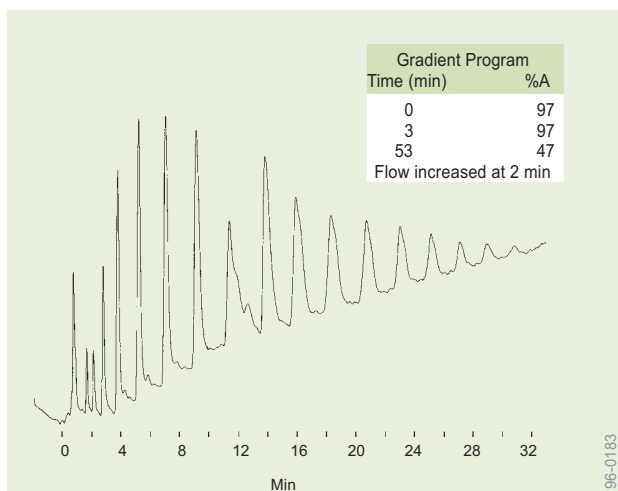

Nonionic Surfactants in Water (SPME/HPLC)

Sample: Triton X-100 in water (500mg/mL)
 SPME Fiber: Carbowax/TPR-100, 65µm film
 Cat. No.: 57315
 Extraction: immersion, 50 min (rapid stirring)
 Desorption: static, 200µL, mobile phase A:B (97:3), 2 min
 Column: SUPELCOSIL LC-NH₂, 25cm x 4.6mm, 5µm particles
 Cat. No.: 58338
 Mobile Phase: A = hexane:2-propanol (90:10) B = 2-propanol:water (90:10)
 Flow Rate: 0.5mL/min (0-2 min) 1.5mL/min (2-53min)
 Det.: UV, 220nm
 Inj.: 200µL

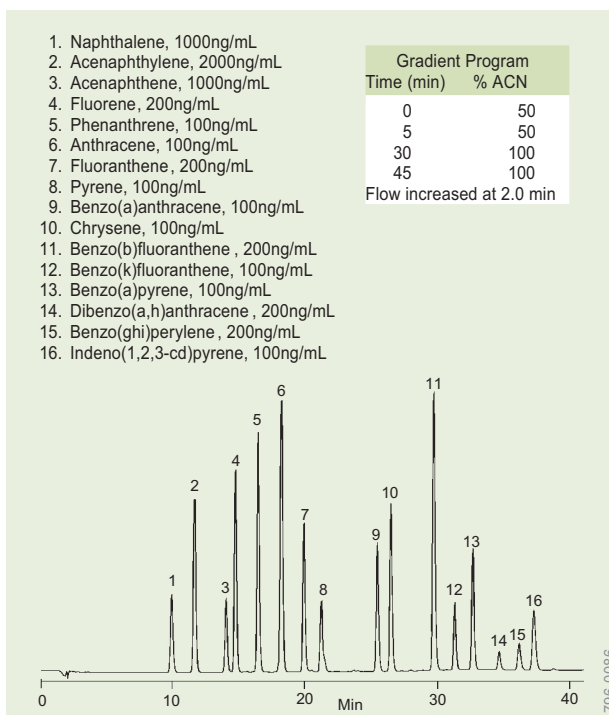

PAHs in Water (SPME/HPLC)

Sample: 5µL PAH mix (Cat. No. 4-8743) in 5mL water
 SPME Fiber: polydimethylsiloxane, 100µm film
 Cat. No.: 57301
 Extraction: immersion, 30 min (rapid stirring)
 Desorption: static, 200µL acetonitrile:water, 40:60, 2 min
 Column: SUPELCOSIL LC-PAH, 15cm x 4.6mm ID, 5µm particles
 Cat. No.: 58318
 Mobile Phase: acetonitrile:water gradient (see program)
 Flow Rate: 0-2 min: 0.2mL/min 2-45 min: 1.0mL/min
 Det.: UV, 254nm


Solid Phase
Microextraction

Order: 1.800.325.3010 Technical Service: 1.800.359.3041 Web: www.sigma-aldrich.com/supelco

SUPELCO

SPME Applications

Semivolatiles, Solvents

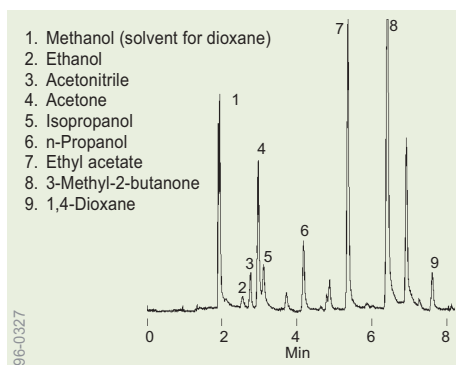

Solvents in Water at 20ppb (SPME /GC)

Sample: solvents at 20ppb in 4mL water + 25% NaCl in 4mL vial

SPME Fiber: Carboxen/PDMS, 75µm film

Cat. No.: 57318

Extractions: immersion, 10 min., rapid stirring

Desorption: 3 min., 260°C

Column: SPB-SULFUR, 30m x 0.32mm ID, 4.0µm film


Cat. No.: 24158

Oven: 50°C (2 min) to 150°C at 10°C/min

Carrier: helium, 30cm/sec

Det.: FID

Inj.: splitless (closed 2 min), 260°C, 0.75mm ID liner

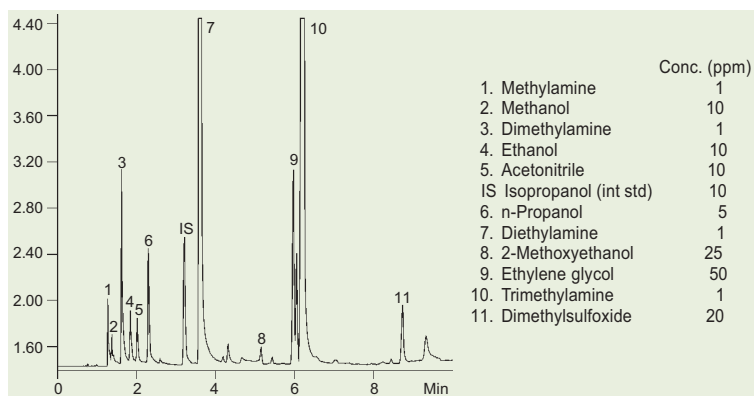

Pharmaceutical Solvents (SPME/GC)

Sample: water + 25% NaCl, pH 11.0 immersion: 3.9mL sample + 0.1mL 50% Na₃PO₄ / 1g NaCl headspace: 2.4mL sample + 0.06mL 50% Na₃PO₄ / 0.6g NaCl

SPME Fiber: polydimethylsiloxane/DVB, 65µm film

Cat. No.: 57310-U

Extraction: immersion, 15 min (rapid stirring or headspace, 15 min, 55°C)

Desorption: 270°C, 5 min

Column: SPB-1 SULFUR, 30m x 0.32mm ID, 4.0µm film


Cat. No.: 24158

Oven: 50°C (2 min) to 180°C at 10°C/min

Carrier: helium, 50cm/sec

Det.: FID

Inj.: splitless (3 min), 270°C (0.75mm ID liner)

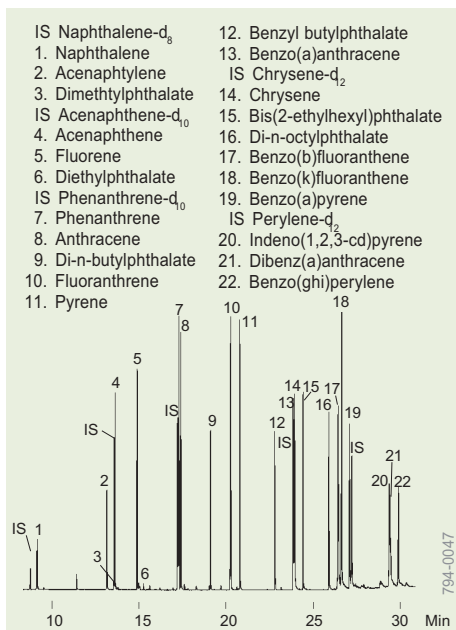

Semivolatiles in Water (SPME/GC)

Sample: water spiked with PAHs and phthalates at 50ppb

SPME Fiber: polydimethylsiloxane, 7µm film

Cat. No.: 57302

Sampling: direct immersion in water, 15 min (rapid stirring)

Inj.: split/splitless, 280°C (closed 4 min)

Column: PTE-5, 30m x 0.25mm ID, 0.25µm film

Cat. No.: 24135-U

Oven: 60°C (3 min) to 320°C at 10°C/min

Carrier: helium, 40cm/sec at 60°C

Det.: MS, Scan Range m/z = 45-465 at 0.6 sec/scan

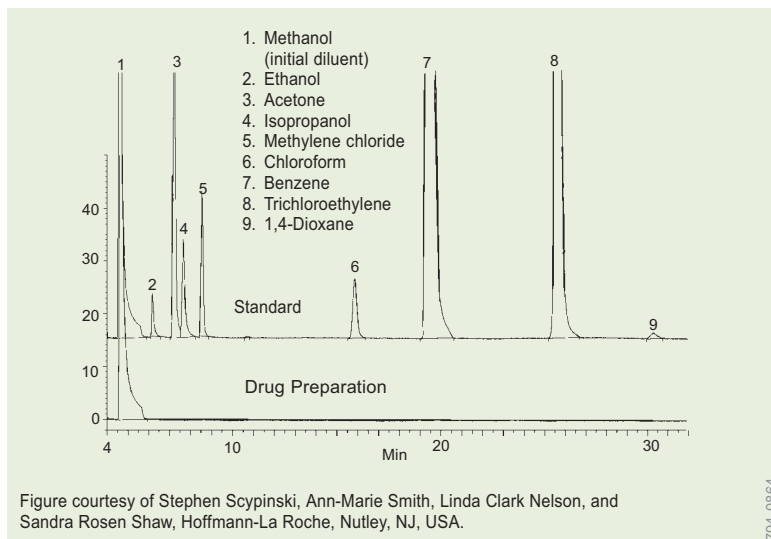


Figure courtesy of Stephen Scypinski, Ann-Marie Smith, Linda Clark Nelson, and Sandra Rosen Shaw, Hoffmann-La Roche, Nutley, NJ, USA.

Residual Solvents in Pharmaceutical Preparations (SPME/GC)

SPME Fiber: polydimethylsiloxane, 100µm film

Cat. No.: 57300-U

Extraction: headspace or immersion sampling (15 min) 1 min desorption

Column: SPB-624, 75m x 0.53mm ID, 3.0µm film

Cat. No.: 25432


Oven: 40°C (35 min) to 220°C at 40°C/min, hold 5 min

Carrier: helium, 35cm/sec

Det.: FID, 250°C


Inj.: splitless (closed 3 min), 200°C

SPME Applications Odors, Sulfur Compounds and Volatiles

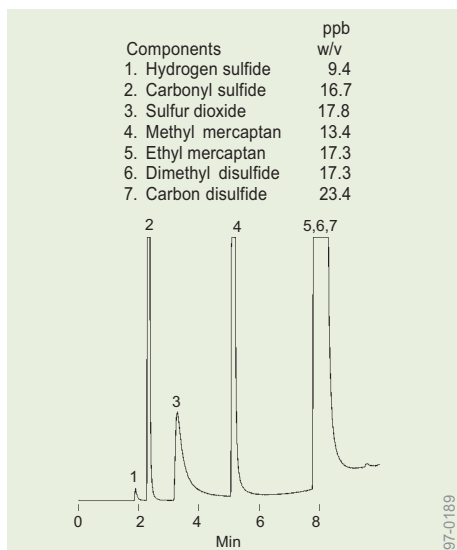


Sulfur Gases (SPME/GC)

Sample: sulfur gases in 40mL vial
 SPME Fiber: Carboxen/PDMS, 75µm film
 Cat. No.: 57318
 Extraction: 10 min headspace at 40°C
 Desorption: 3 min at 290°C
 Column: Supel-Q PLOT, 30m x 0.32mm ID
 Cat. No.: 24242
 Oven: 45°C (1 min) to 250°C at 15°C/min

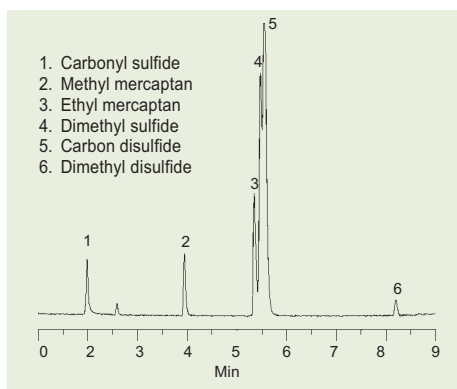


Sulfur Gases at 1ppm (SPME/GC)

Sample: sulfur gases in air at 1ppm, 250mL bulb
 SPME Fiber: Carboxen/PDMS, 75µm film
 Cat. No.: 57318
 Extraction: headspace, 5 min., ambient temp.
 Desorption: 2 min., 250°C
 Column: Supel-Q PLOT, 30m x 0.32mm ID
 Cat. No.: 24242
 Oven: 45°C (0.75 min.) to 250°C at 25°C/min
 Carrier: helium, 25cm/sec
 Injection: splitless/split (closed 2 min.), 0.75mm ID liner
 Detector: MS quadrupole, m/z = 32 - 125 (0.6sec/scan)

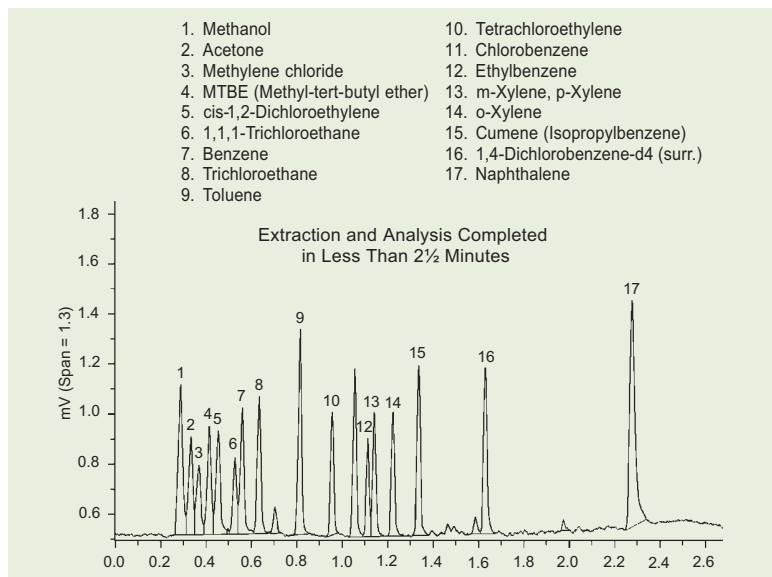

Water Sample Screened for BTEX and Other Compounds (SPME/GC)

Sample: 0.7mL water spiked with 2-600µg/liter each analyte, 0.25g NaCl added
 SPME Fiber: 30µm PDMS
 Cat. No.: 57309
 Extraction: headspace, 12 sec (no stirring), ambient temp.
 Desorption: 2-3 min, 250°C (0.75-1.0mm ID inj. port liner)
 GC Column: Equity-1, 10m x 0.20mm ID, 1.2µm film (available on request)
 Oven: 70°C (0.2 min) to 180°C at 50°C/min
 Carrier: hydrogen, 12psi (+ 30mL/min nitrogen make-up gas)
 Det.: FID, 250°C

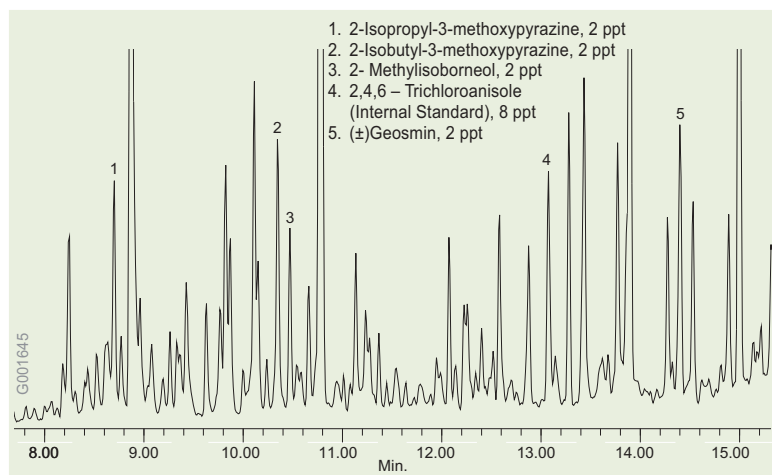

Trace Odors in Drinking Water (SPME/GC)


Sample: 25mL of water containing 25% NaCl and drinking water odors kit, Cat. No. 47529-U
 SPME Fiber: 2cm StableFlex coated with 50/30µm DVB/Carboxen/PDMS
 Cat. No.: 57348-U
 Extraction: headspace, 65°C (30 min.)
 Column: Equity-5, 30m x 0.25mm, 0.25µm film
 Cat. No.: 28089-U
 Desorption: 3 min. at 260°C
 Oven: 60°C (2min) to 200°C at 8°C/min
 GC Liner: 0.75mm SPME liner
 Detector: MS, selected ions (SIM) 95, 112, 124, 137, 197; interface at 280°C
 Flow: Helium, 37cm/sec @ 60°C (1mL/min constant flow)
 Inj.: SPME fiber, splitless opened at after 1 min at 50mL/min.

SPME Applications

Volatiles

Solid Phase
Microextraction

Order: 1.800.325.3010 Technical Service: 1.800.359.3041 Web: www.sigma-aldrich.com/supelco

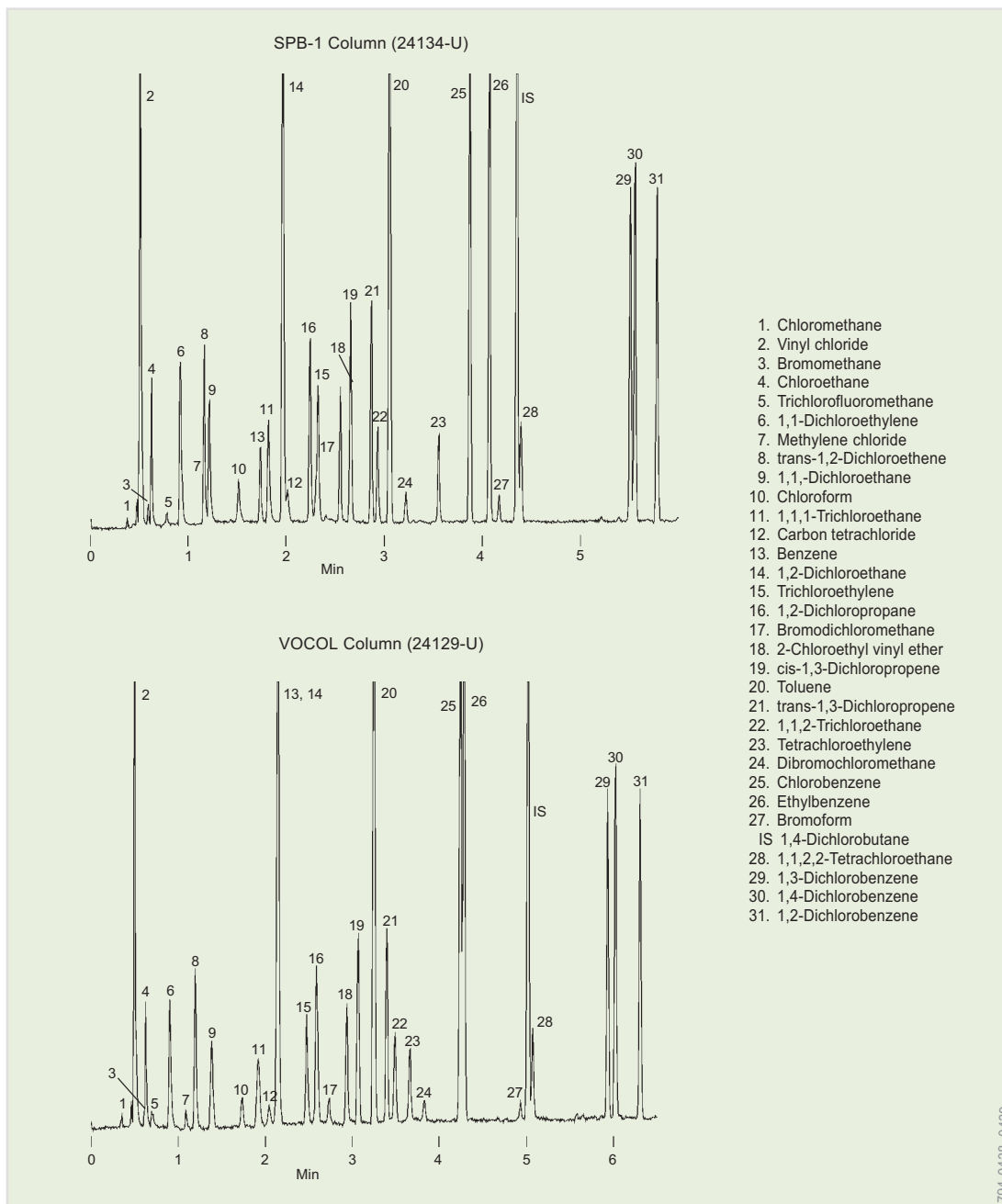


794-0438, 0439

Volatiles (SPME/GC)

Sample: 50ppb each analyte in 1.8mL saturated salt water
 SPME Fiber: polydimethylsiloxane, 100µm film
 Cat. No.: 57300-U
 Extraction: 5 min immersed in water with rapid stirring
 Columns: 10m x 0.20mm ID, 1.2µm film
 Oven: 40°C (0.75 min) to 160°C at 20°C/min
 Carrier: helium, 40cm/sec
 Det.: FID, 260°C
 Inj.: 230°C, splitless (closed 3 min)

- | | | | |
|-------------------------------|-------------------------------|-------------------------------|---------------------------------------|
| 1. Dichlorodifluoromethane | 17. Carbon tetrachloride | 32. Chlorobenzene | 47. 4-Chlorotoluene |
| 2. Chloromethane | 18. 1,2-Dichloroethane | 33. 1,1,1,2-Tetrachloroethane | 48. tert-Butylbenzene |
| 3. Vinyl chloride | 19. Benzene | 34. Ethylbenzene | 49. 1,2,4-Trimethylbenzene |
| 4. Bromomethane | IS Fluorobenzene | 35. m-Xylene | 50. sec-Butylbenzene |
| 5. Chloroethane | 20. Trichloroethylene | 36. p-Xylene | 51. p-Isopropyltoluene |
| 6. Trichlorofluoromethane | 21. 1,2-Dichloropropane | 37. o-Xylene | 52. 1,3-Dichlorobenzene |
| 7. 1,1-Dichloroethylene | 22. Bromodichloromethane | 38. Styrene | 53. 1,4-Dichlorobenzene |
| 8. Methylene chloride | 23. Dibromomethane | 39. Isopropylbenzene | 54. n-Butylbenzene |
| 9. trans-1,2-Dichloroethylene | 24. cis-1,3-Dichloropropene | 40. Bromoform | IS 1,2-Dichlorobenzene-d ₄ |
| 10. 1,1,-Dichloroethane | 25. Toluene | 41. 1,1,2,2-Tetrachloroethane | 55. 1,2-Dichlorobenzene |
| 11. 2,2-Dichloropropane | 26. trans-1,3-Dichloropropene | 42. 1,2,3-Trichloropropane | 56. 1,2-Dibromo-3-chloropropane |
| 12. cis-1,2-Dichloroethylene | 27. 1,1,2-Trichloroethane | 43. n-Propylbenzene | 57. 1,2,4-Trichlorobenzene |
| 13. Chloroform | 28. 1,3-Dichloropropane | 44. Bromobenzene | 58. Hexachlorobutadiene |
| 14. Bromochloromethane | 29. Tetrachloroethylene | 45. 1,3,5-Trimethylbenzene | 59. Naphthalene |
| 15. 1,1,1-Trichloroethane | 30. Chlorodibromomethane | 46. 2-Chlorotoluene | 60. 1,2,3-Trichlorobenzene |
| 16. 1,1-Dichloropropene | 31. 1,2-Dibromoethane | | |

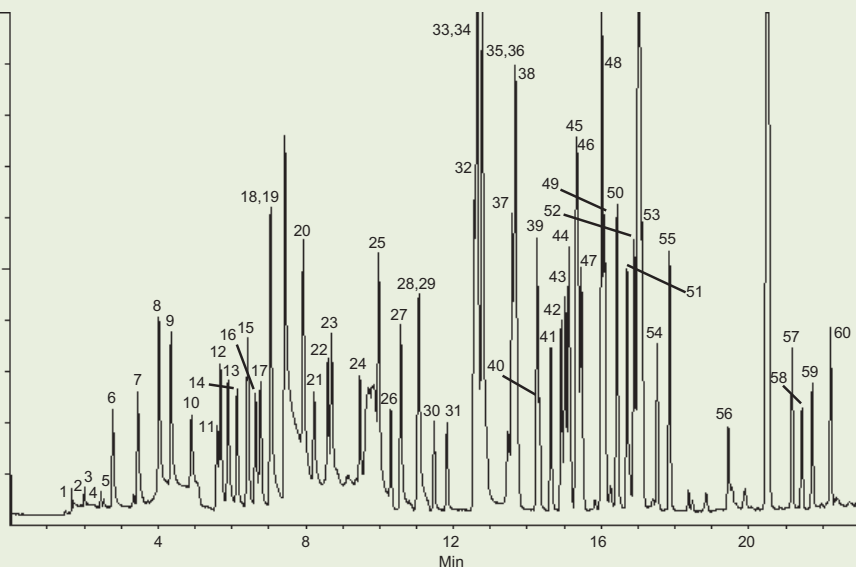

97-0183

Volatile Compounds by US EPA 524.2 (SPME/GC)

Sample: 1ppb each analyte in 3mL water (25% NaCl added) in 4mL vial
 SPME Fiber: Carboxen/PDMS, 75µm film
 Cat. No.: 57318
 Extraction: headspace, 40°C (20 min.), with stirring
 Desorption: 5 min., 310°C
 Column: VOCOL, 30m x 0.25mm ID, 1.5µm film
 Cat. No.: 24205-U
 Oven: 40°C (2 min.) to 210°C at 8°C/min
 Carrier: helium, 35cm/sec
 Injection: splitless/split (closed 2 min), 0.75mm ID liner
 Detector: ion trap, MS, m/z = 45 - 260 (0.6sec/scan)